

For Windows10, please note the followings:

When installing coolingtech measurement, please right click the install file Measurement.exe, select 'run as administrator'.

when using coolingtech measurement software, please do as the followings:

- 1, After installing the software, find "CoolingTech_Microscope" by the installation file folder path ,if you choose the default path, please select the file folder "CoolingTech_Microscope" under the file folder "C:\Program Files (x86)\CoolingTech";
- 2, Right click on the folder "CoolingTech_Microscope" and select "properties" - > "safe" - > "Edit" button, and then click "add (D)" button, please enter "everyone" under the blank of "enter the object name to choose:" , and then click "OK" button;
- 3 , Please click "Everyone" under "group or user name (G):" , --> "Permission of Everyone" --> "complete control" --> "allow";
- 4, click on the "application" and "OK" button .

then the coolingtechmeasurement software can be used normally.